

# Algae Online Monitor AOM 2700

# Algae Online Monitor AOM 2800

## Operation Manual



Please read this guide before operating the AOM device

Photon Systems Instruments, Högrova 20, 612 00 Brno,  
Czech Republic, <http://www.psi.cz>

# Table of Contents

---

---

- 1. List of Equipment..... 3**
- 2. General Information ..... 4**
- 3. Technical Specification ..... 5**
- 4. Maintenance and General Safety Guidelines..... 7**
- 5. Physical Features..... 8**
- 6. Cable Connections..... 9**
- 7. Device Installation ..... 11**
- 8. Operation Instructions..... 13**
- 9. FluorPen Software..... 21**
  - 9.A. Menu and Icon Explanation..... 22*
  - 9.B. Example of Data Transfer and Visualization..... 24*
  - 9.C. Explanation of OJIP Parameters..... 26*
  - 9.D. Software Update ..... 27*
- 10. Statement of Limited Warranty..... 29**

# 1. List of Equipment

---

Carefully unpack the carton. You should have received the following items:

- **Algae Online Monitor 2700 or Algae Online Monitor 2800**
- **Flow-Through Cuvette**
- **Anti-Vibration Pad**
- **Mounting Clips**
- **Cables and Connectors**
- **This Operation Manual** (*on a CD or a printed version*)
- **Install CD with the FluorPen software**
- **Optional Accessories (Hoses, Supplementary Cuvette, Pump, Control Valve**  
*(according to your specific order)*

Note:

If any item is missing, please, contact the manufacturer. Also check the carton for any visible external damage. If you find any damage, notify the carrier and the manufacturer immediately. The carton and all packing materials should be retained for inspection by the carrier or insurer.

For customer support, please write to: [support@psi.cz](mailto:support@psi.cz)

## 2. General Information

---

Algae Online Monitor (AOM) is a portable, robust device for online detection and continuous monitoring of photosynthetic microorganisms in both natural and artificial water bodies. It can serve for detection and discrimination among variety of cyanobacteria, green and brown algae, diatoms, and other microbes. Its ultra-high sensitivity (30 ng Chl/l) allows early detection of very low concentrations of these organisms.

According to customer's needs, the AOM is equipped with two measuring lights:

- Either with blue (455 nm) and red (630 nm) light.
  - Or with blue (455 nm) and amber (590 nm) light.
- 

**The AOM device measures the following chlorophyll fluorescence parameters:**

### **$F_T$ - Instantaneous Fluorescence**

$F_T$  is equivalent to  $F_0$  if the sample is dark-adapted.  $F_T$  is equivalent to  $F_S$  in a light-adapted sample.

### **QY - Quantum Yield**

QY is a measure of the Photosystem II efficiency. In a dark-adapted sample this is equivalent to  $F_v/F_m$ . In a light-adapted sample it is equivalent to  $F_v'/F_m'$ .

### **OJIP - Chlorophyll Fluorescence Induction Kinetics**

OJIP transients enable observing major changes that occur during exposure of a sample to high irradiance. This biophysical signal is extremely rich in terms of information and reflects the time course of photosynthesis. In AOM device, OJIP is represented by the Fix Area Parameter.

Fix Area = Total area above the OJIP fluorescence transient. This parameter correlates with total change of fluorescence signal in OJIP protocol, with total pigment content and hence with cell concentration.

### 3. Technical Specification

---

---

#### Measured and Calculated Parameters:

$F_T$ , QY, OJIP (Fix Area)

#### Detection Limit:

Algae – 10 cells/ml; cyanobacteria – 100 cells/ml

#### Detector Wavelength Range:

PIN photodiode with 667 to 750 nm bandpass filters

#### Actinic and Saturating Light:

Adjustable from 0 to 3,000  $\mu\text{mol}(\text{photons})/\text{m}^2.\text{s}$

#### Measuring Light:

Blue (455 nm) and red (630 nm) measuring light adjustable by intensity

Blue (455 nm) and amber (590 nm) measuring light adjustable by intensity

#### Communication:

Serial RS 232, RS 485

#### FluorPen Software:

Windows 2000, XP, or higher\*

#### Bios:

Upgradeable firmware

#### Memory Capacity:

4 Mb – up to 100,000 data points (about 300 OJIP curves)

#### Display:

AOM 2800: 2 x 16 characters LC display

AOM 2700: without display

#### Keypad:

AOM 2800: 2-key tactile response

AOM 2700: remote control

#### Power Saving Mode:

Automatic

#### Power Supply:

Power source 24 V (optionally 12 V)

#### Size:

20 cm x 23 cm x 11 cm

#### Weight:

3.4 kg

#### Sample Holder:

Flow-through cuvette made of quartz glass

\* Windows is a registered trademark of Microsoft Corporation.

**Case:**

**Waterproof IP65**

**Operating Conditions:**

**Temperature: 0 to 45 °C; Relative humidity: 0 to 95 % (non-condensing)**

**Warranty:**

**1 year parts and labor (see the last page of this Operation Manual for precise warranty conditions)**

## 4. Maintenance and General Safety Guidelines

---

- **Connect all peripheral units when the device is switched off!**
- **First connect all peripheral units (pumps, control valves), then switch on the device.**
- **Never submerge the whole device in water!**
- Keep the optical part clean and dry. If cleaning is needed, use the soapy water and soft, non-abrasive tissue.
- The device should not come in contact with any organic solvents, strong acids or bases.


### **GENERAL ELECTRICAL SAFETY GUIDELINES:**

- Routinely check the devices and their wiring.
- Replace worn or damaged cords immediately.
- Use electrical extension cords wisely and do not overload them.
- Place the devices on a flat and firm surface. Keep them away from wet floors and counters.
- Avoid touching the device, socket outlet or switch if your hands are wet.
- Do not perform any alterations to the electrical part of the devices or its components.

## 5. Physical Features

---


### Algae Online Monitor AOM 2700

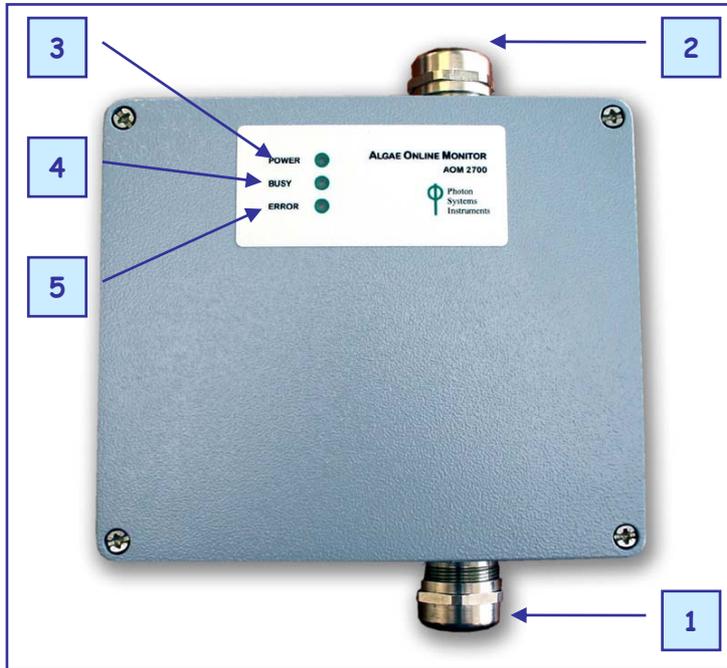

- [1] Inlet
- [2] Outlet

Three LED Indicators:  
[3] On / Off  
[4] Device operation  
[5] Error

### Algae Online Monitor AOM 2800

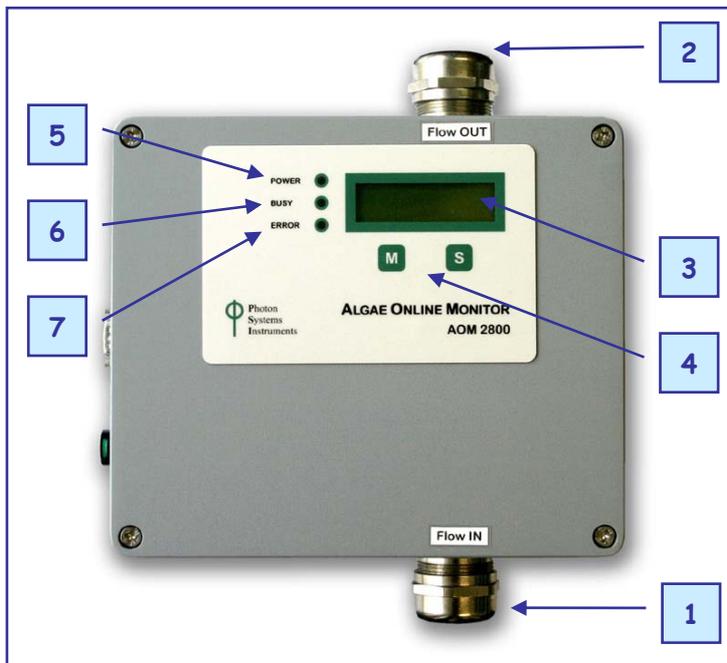


- [1] Inlet
- [2] Outlet
- [3] Display
- [4] Control keys

Three LED Indicators:  
[5] On / Off  
[6] Device operation  
[7] Error

## 6. Cable Connections

---

### Algae Online Monitor AOM 2700

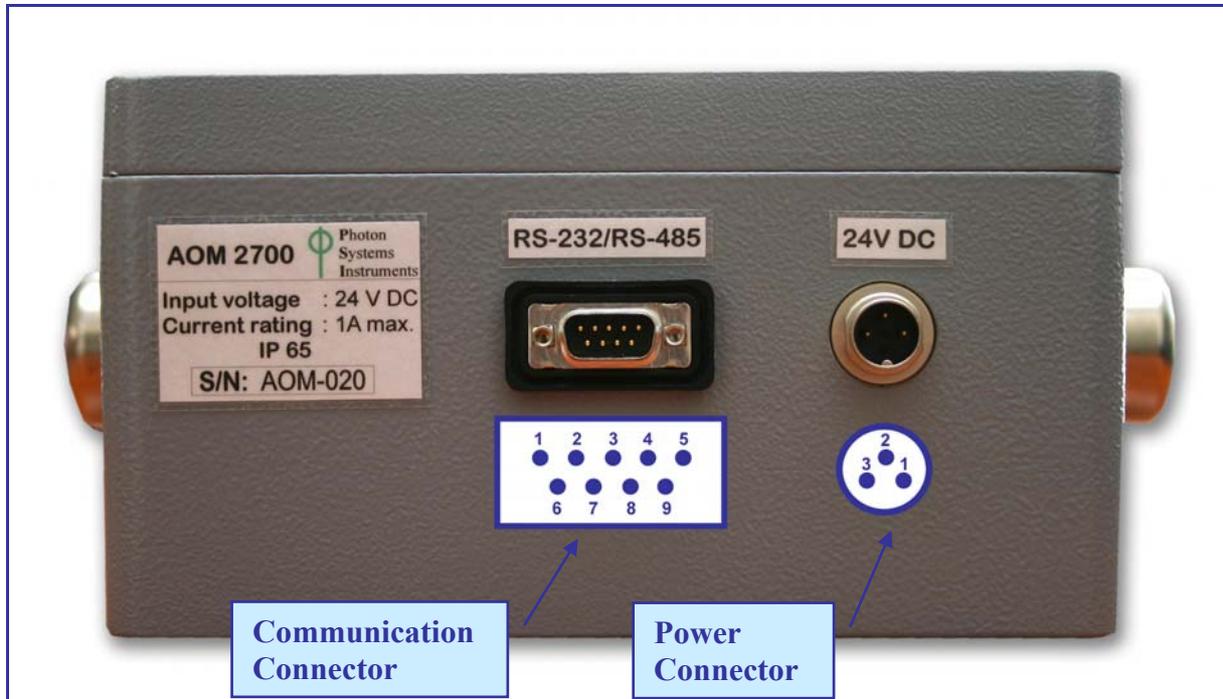

#### Communication Connector:


Serial RS-232: [2] - RXD  
                  [3] - TXD  
                  [5] - GND

Serial RS-485: [6] - Y  
                  [7] - Z  
                  [8] - A  
                  [9] - B

#### Power Connector:

[1] - +24 V, +12 V  
[2] - GND

## Algae Online Monitor AOM 2800

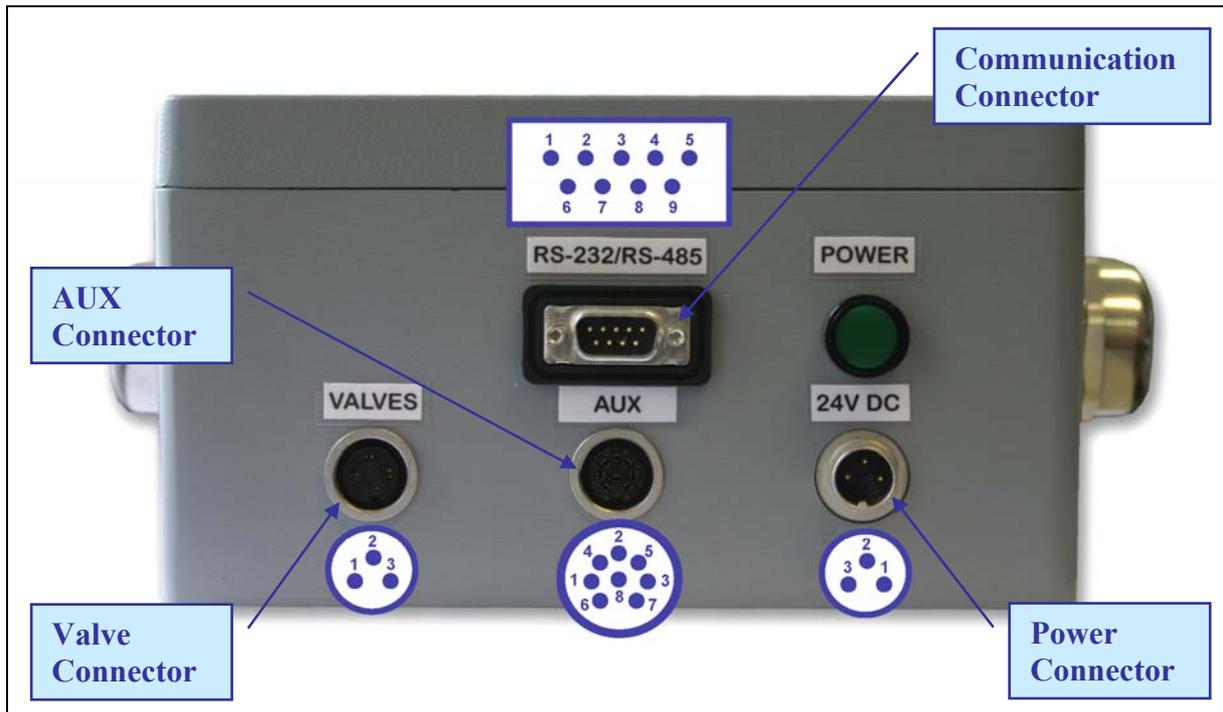

### Communication Connector:

Serial RS-232: [2] - RXD  
[3] - TXD  
[5] - GND

Serial RS-485: [6] - Y  
[7] - Z  
[8] - A  
[9] - B

### Power Connector:

[1] - +24 V, +12 V  
[2] - GND

### Valve Connector:

[1] - +24 V, 12 V, Common Valve  
[2] - Main Valve/Pump  
[3] - Cleaning Valve/Pump

### AUX Connector:


[1] - Current Loop  
[2] - GND  
[3], [4], [5], [6] - Digital Out  
[7] - Alarm  
[8] - +24V, +12 V

## 7. Device Installation

---


### 1. Installation of the measuring instrument on a wall (or other base)

- A) fix four elastic elements(1) on the wall
- B) use four stainless steel nuts(2) and fix the device

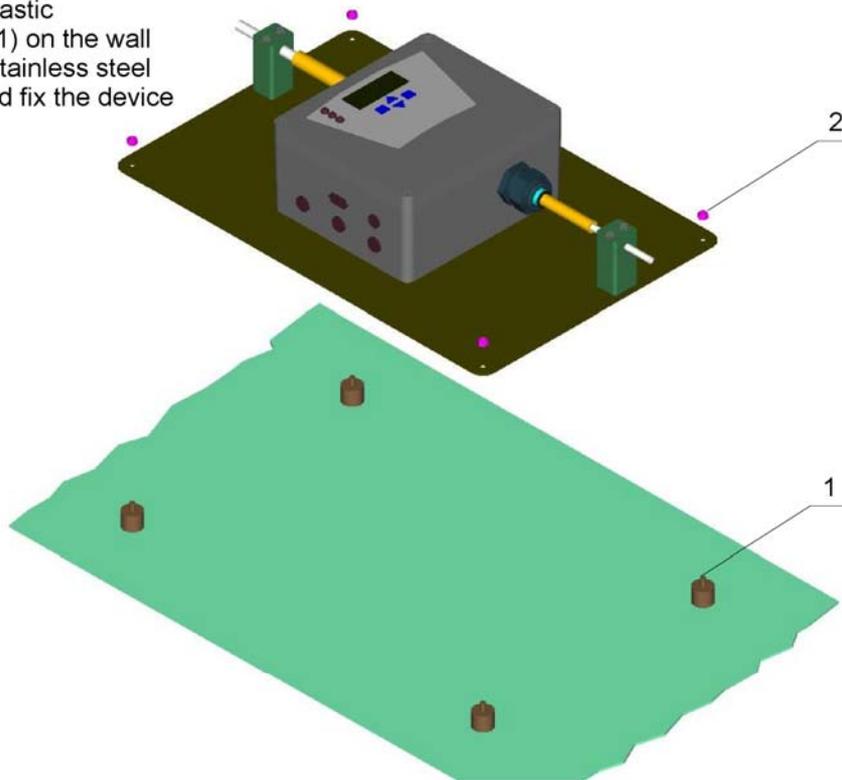


### 2. Optimal configuration of the measuring instrument

- (3) input tube
- (4) output tube

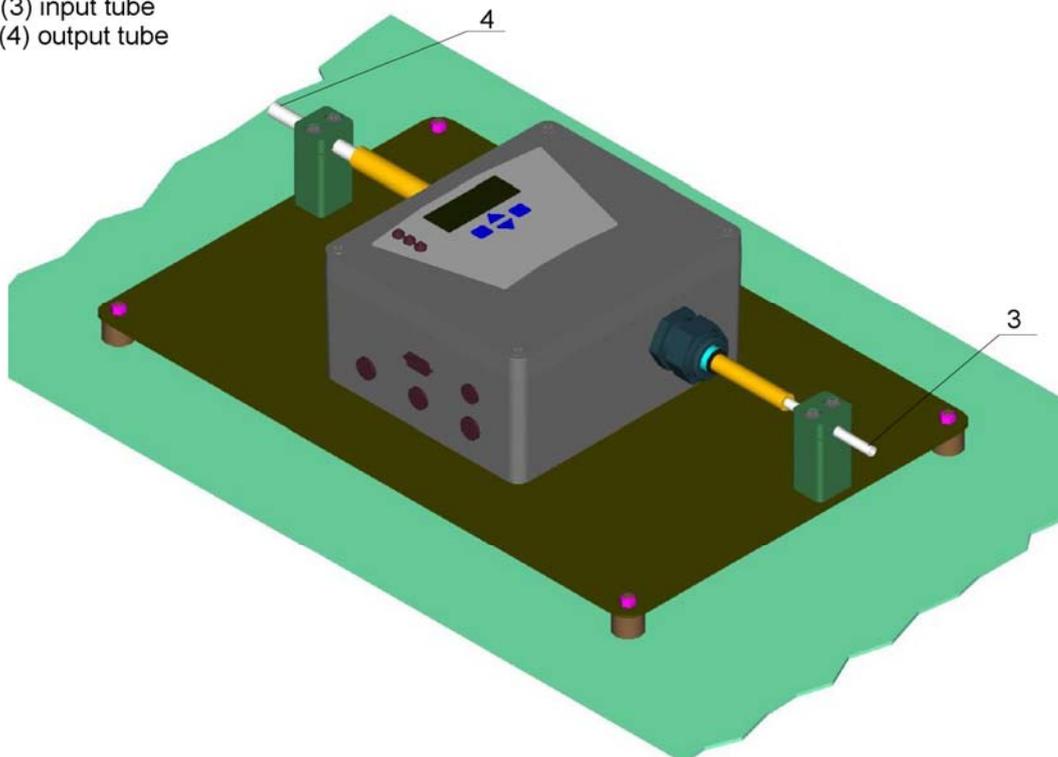


### 3. Installation of the input and output tubes

- (3) input tube
- (4) output tube

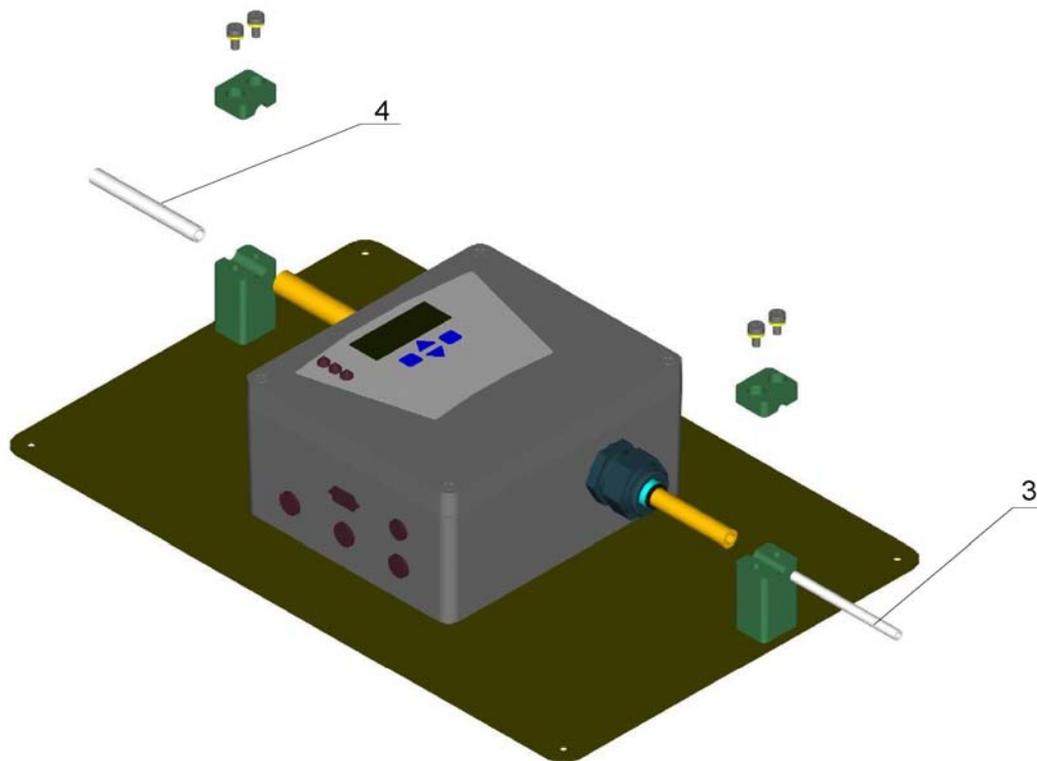

### 4. Replacement of the flow-through cuvette (5)

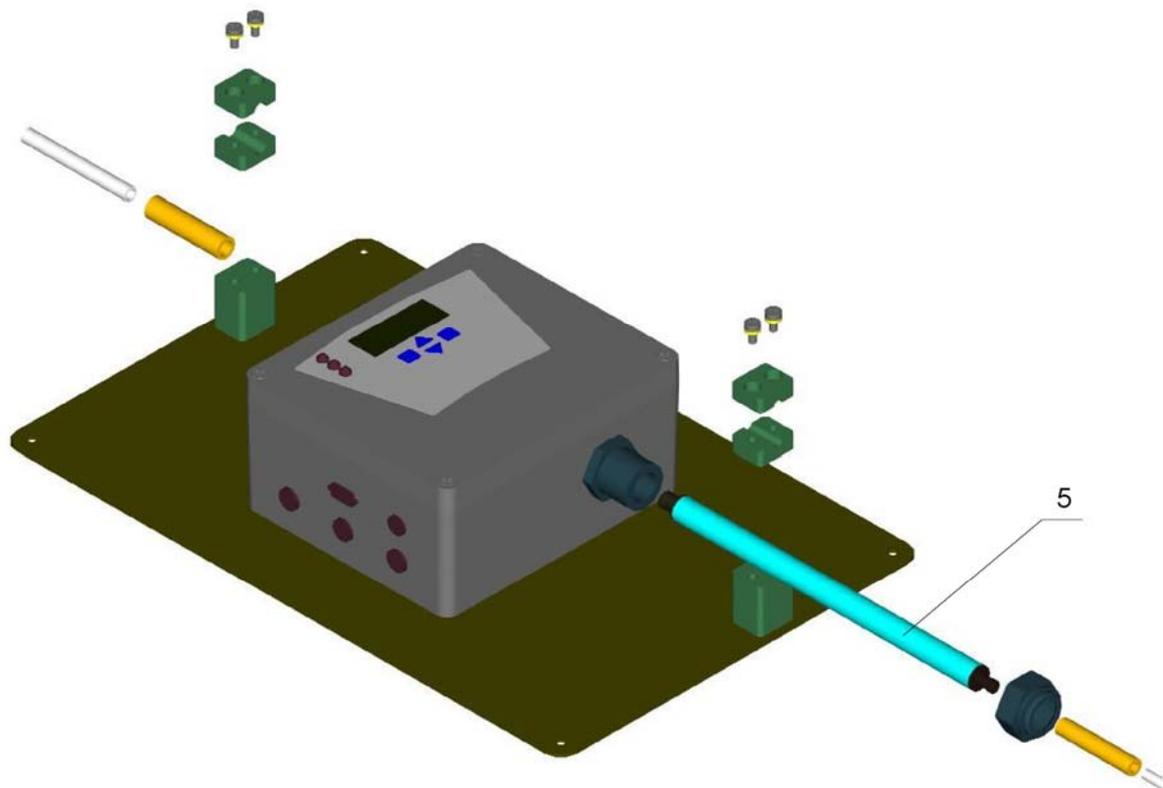

## 8. Operation Instructions

---

**Algae Online Monitor is operated by two keys:**

- “S” key allows the user to select a menu option (based on cursor > position), or to move forward to a “lower level” position.
- “M” key allows the user to scroll through single menu and sub-menu options, or to move back to a “higher level” position.

**The next six pages bring graphical presentation of the operation scheme, which is structured into three levels:**


- **Main Menu**
- **First-level Sub-Menus** (Measure, Cleaning, Data Browse and Erase, Setting, Time Indication)
- **Second-level Sub-Menus** (see next pages)

**Explanation of symbols and color differentiation\* used in the graphical presentation:**


- The **blue** color represents the **Main Menu** and its Options.
- The **yellow** color represents the **first-level Sub-Menus** and their Options.
- The **green** color represents the **second-level Sub-Menus** and their Options.
- Full-line arrows are used for the “S” key. 
- Dashed-line arrows are used for the “M” key. 

*\* The AOM display does not reflect this color differentiation.*

# Main Menu

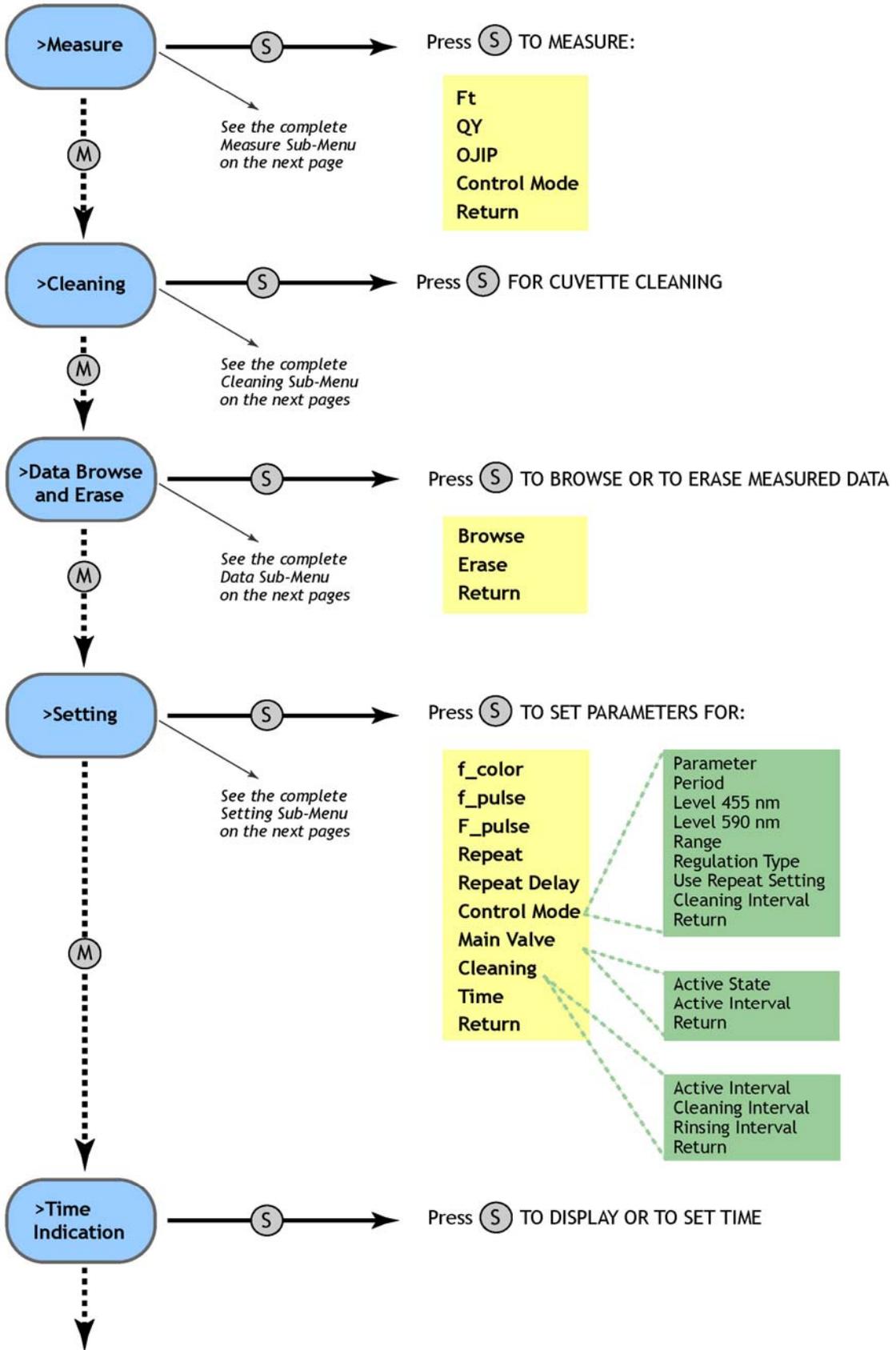


## Measure Sub-Menu

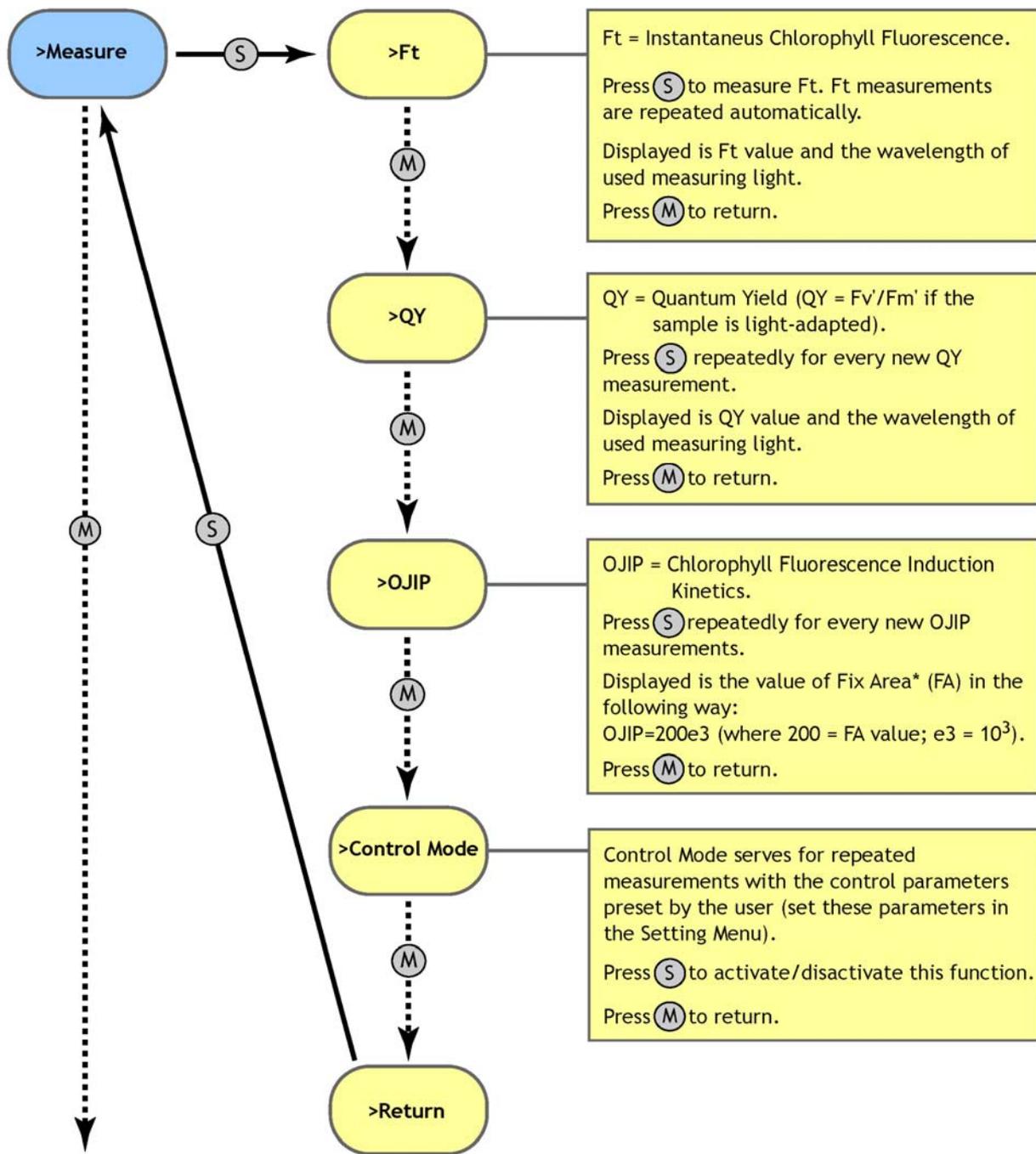


\* Fix Area = Total area above the OJIP fluorescence transient and total change of fluorescence signal in OJIP protocol. This parameter correlates with total pigment content and hence with cell concentration.


## Cleaning and Data Sub-Menu

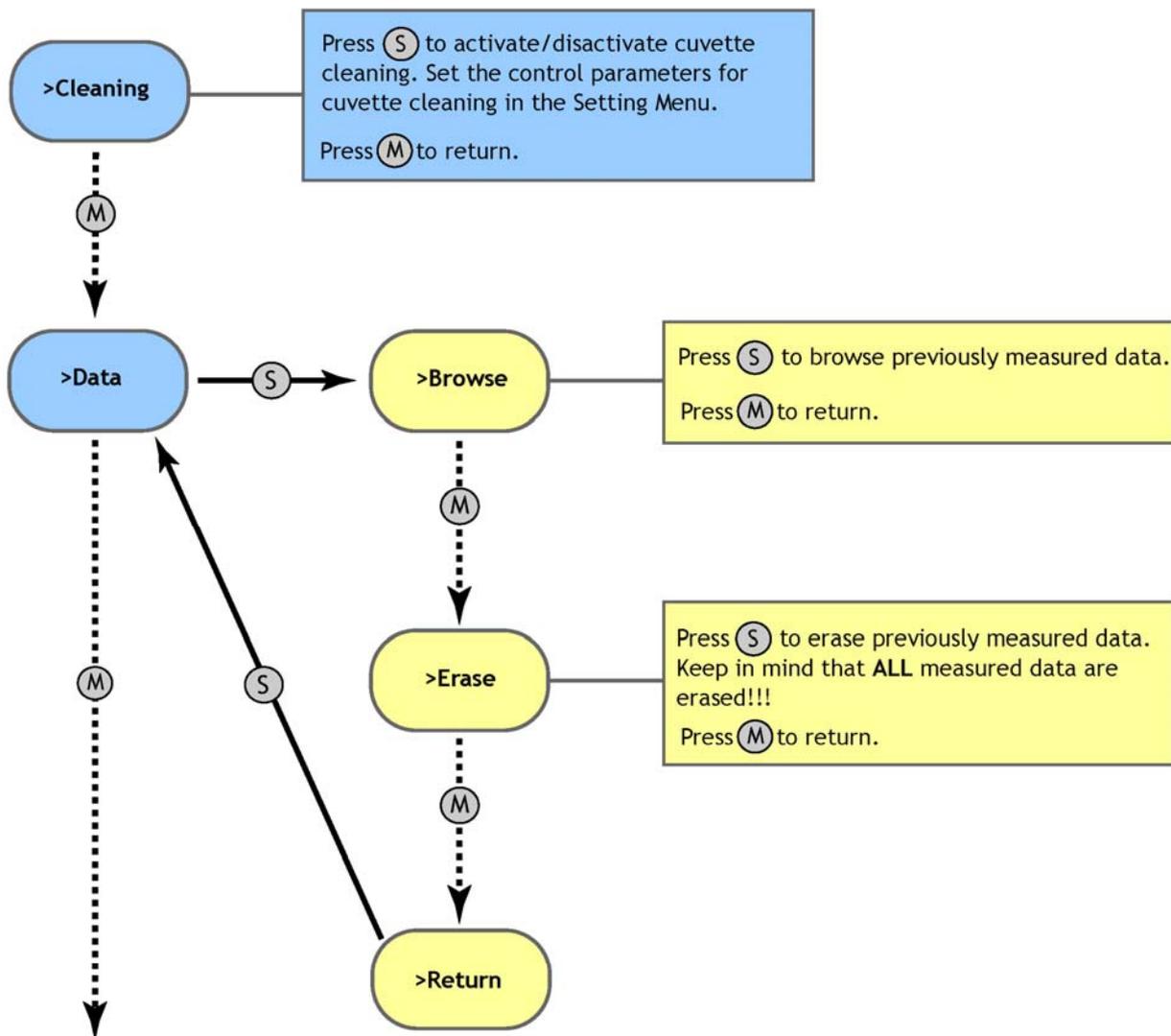


## Setting Sub-Menu (Part 1)

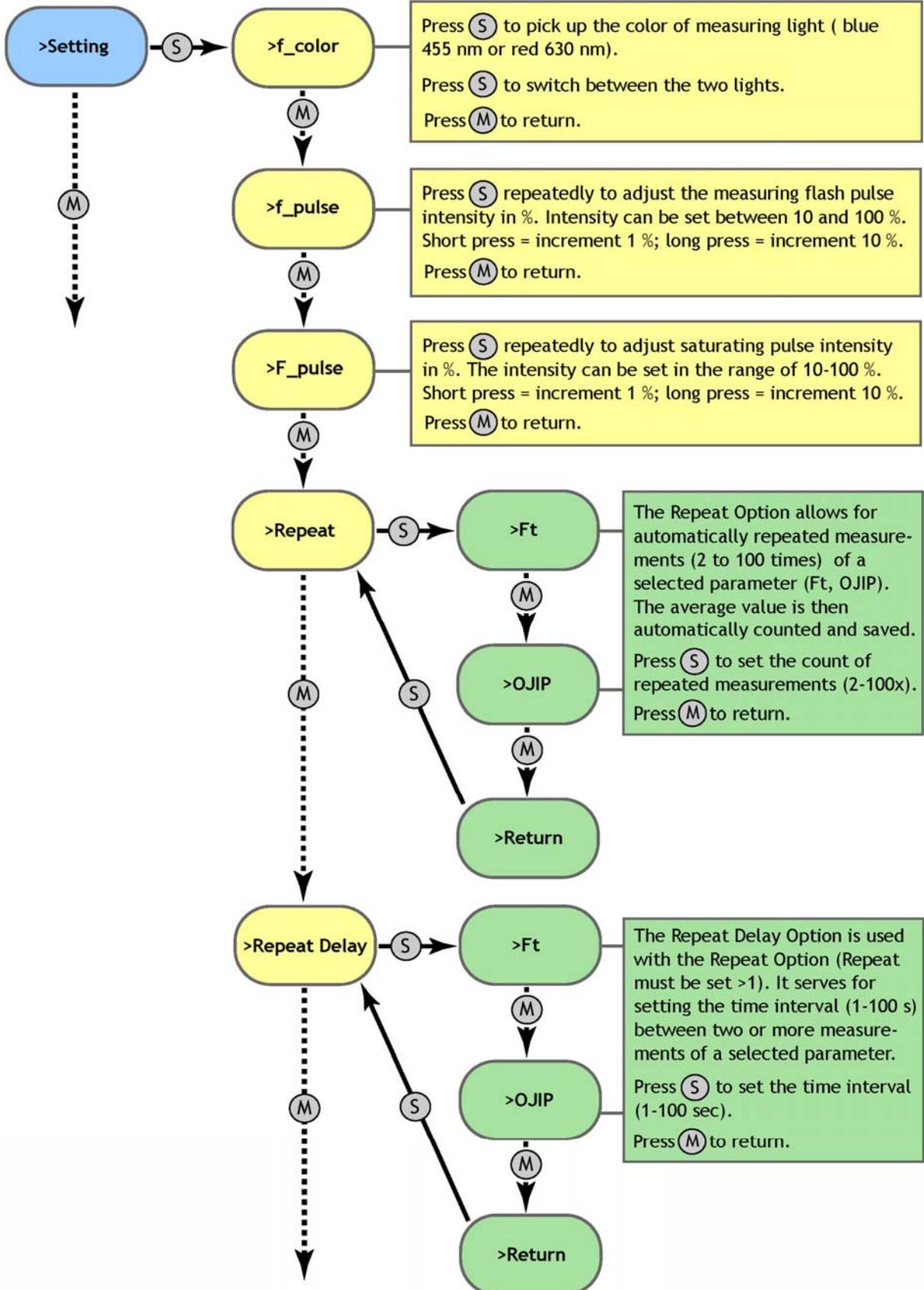


## Setting Sub-Menu (Part 2)

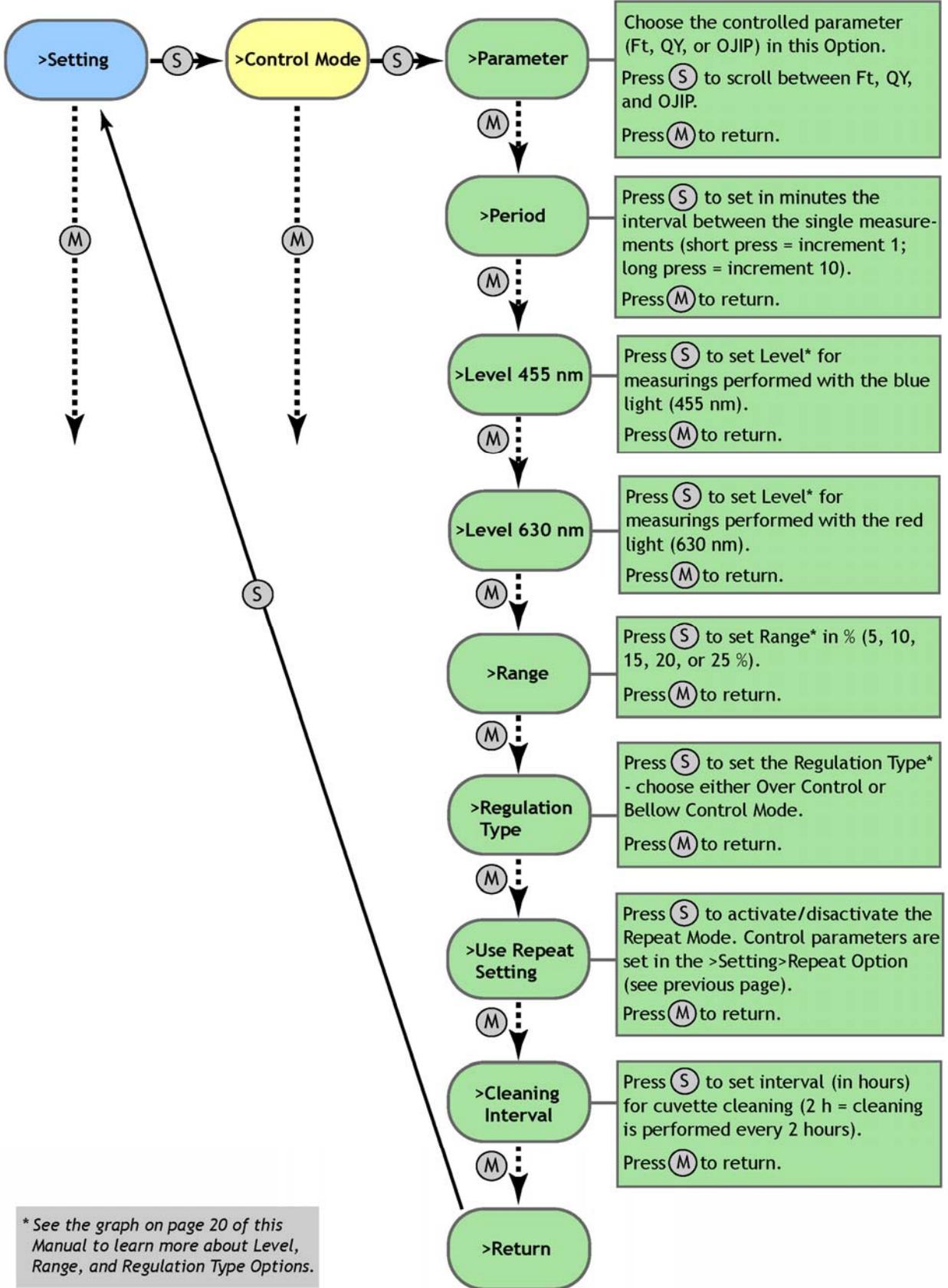

## Setting Sub-Menu (Part 3)

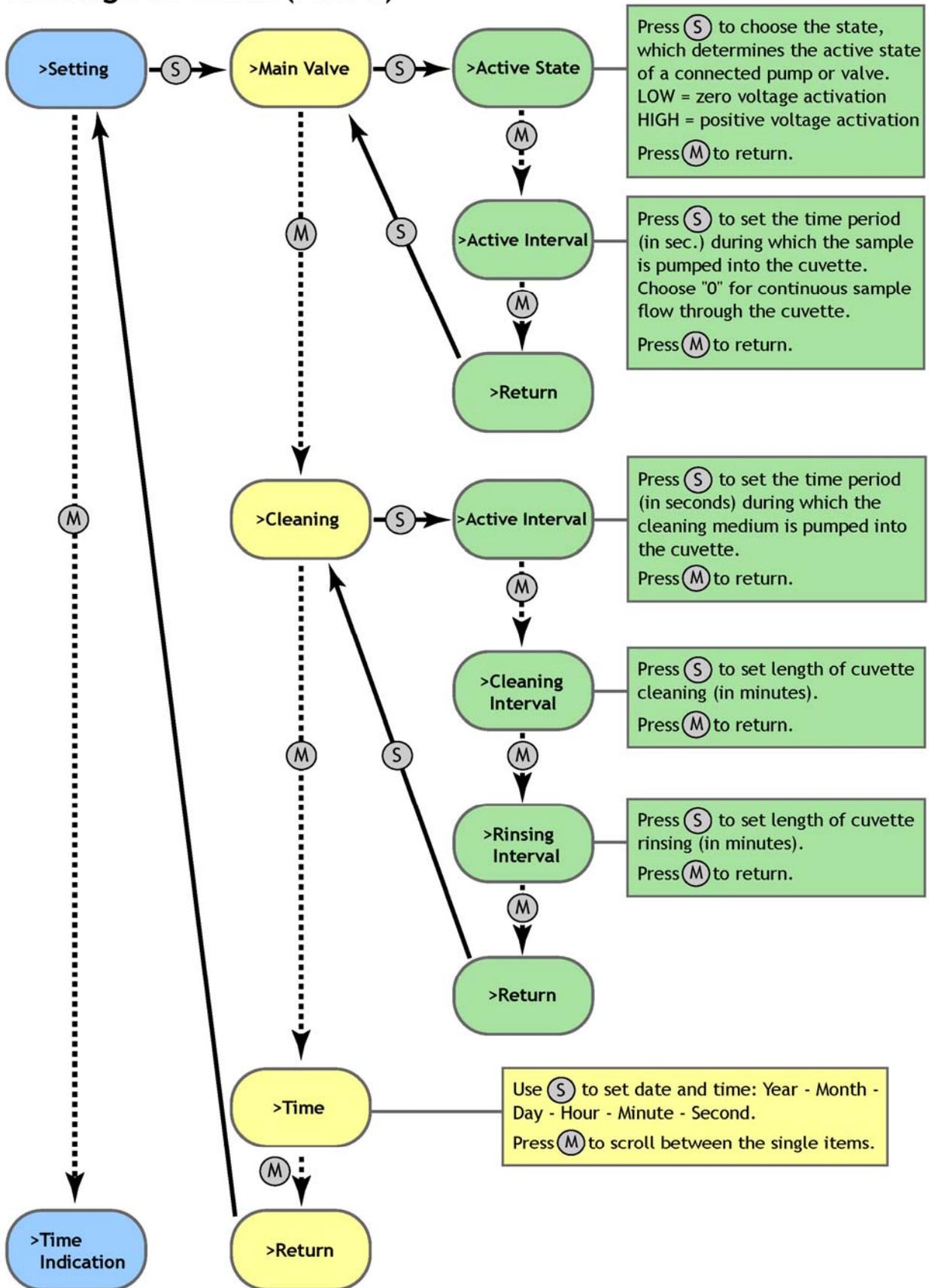

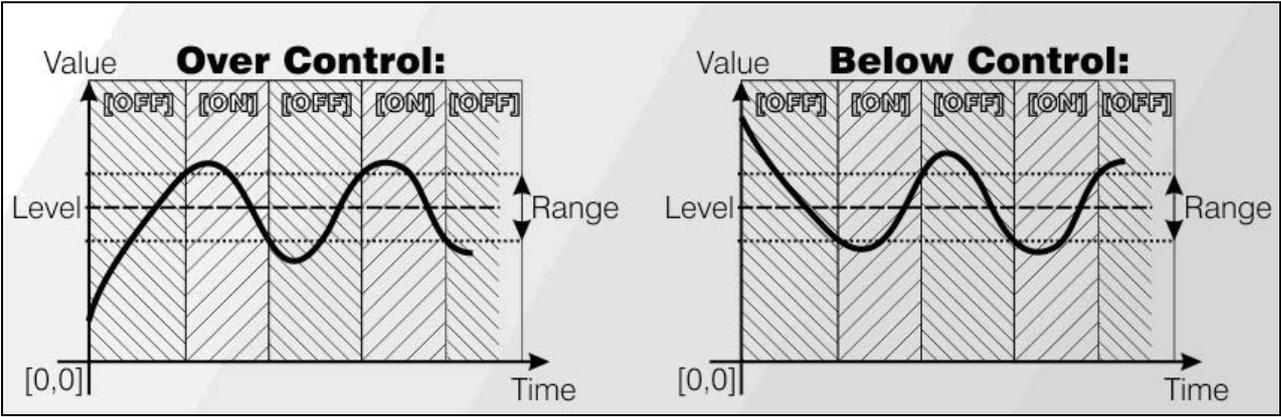

## 9. FluorPen Software

---


### Starting up:

1. Switch on the computer.
2. Switch on the AOM device.
3. Start the FluorPen program.

### Connecting:

**Select:** Setup>Device ID (Ctrl+I).

If properly connected, the message “Device: FluorPen” appears in the bottom part of the screen.

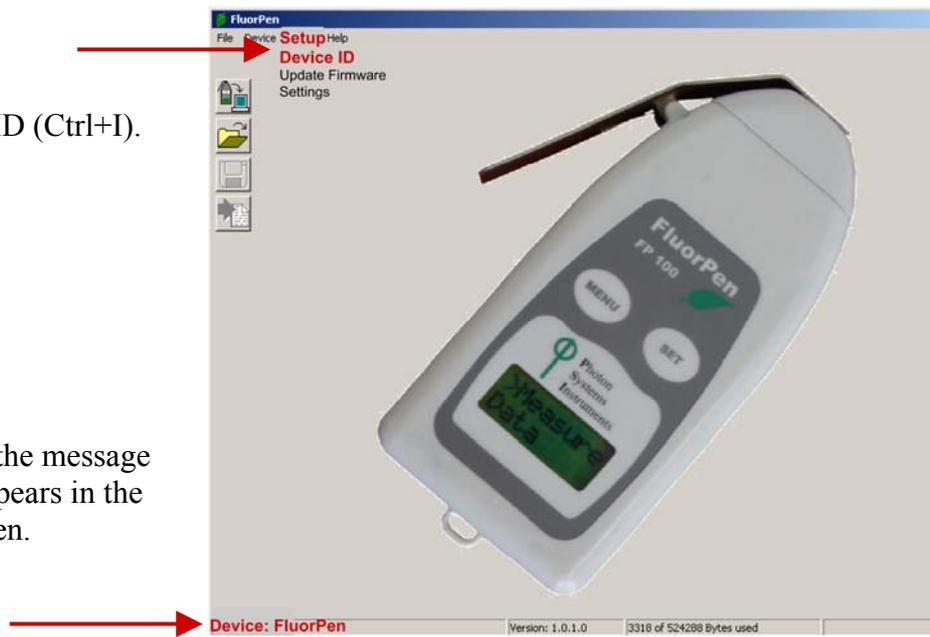


## 9.A. Menu and Icon Explanation

---


### Menu: File

- Load** Loads previously saved data files.
- Save** Saves data to hard disc.
- Export** Exports data in .txt format.
- Close** Closes the current experiment.
- Close All** Closes all running experiments.
- Exit** Exits the program.

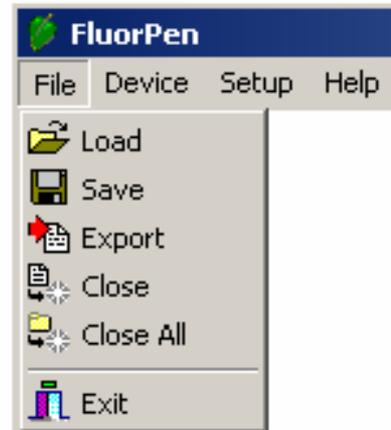

### Menu: Device

- Download** Downloads data from the AOM to your PC.
- Erase Memory** Erases data from the AOM memory.

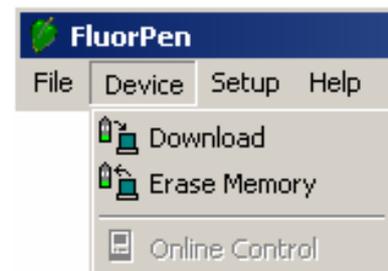

### Menu: Setup

- Device ID** Detects the connected device.
- Update Firmware** Used for software updates.\*
- Settings** Used for modification of the program settings.\*\*




\* For more information on software updating, see Chapter 8.D. of this Operation Manual.

\*\* See more information on the next page.

### Menu: Help

- About** Offers basic information about the program.
- Register** Used for the FluorPen software registration.

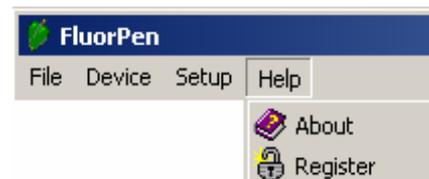

---

## Menu: Settings

### After Download - Memory Erase


If the box is checked the AOM memory is erased after each data download.

### Data - Inverted

If the box is checked the polarity of data is inverted, e.g., multiplied by -1.\*

### Data - Add to Opened

If the box is checked the downloaded data are added to that of the current opened experiment.

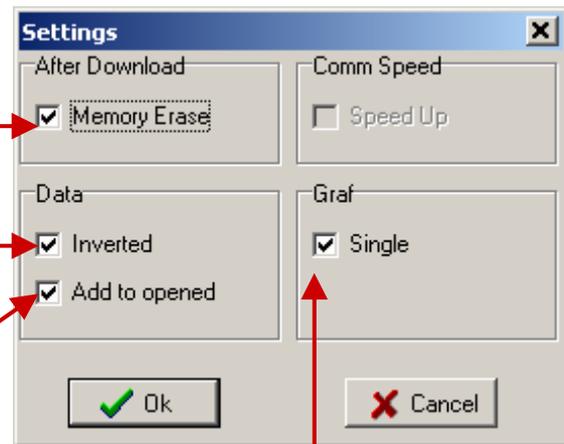

### Graf - Single


If the box is checked all measured data are visualized in one graph, i.e., the value of each new measurement is added to the currently used graph window.

If the box is not checked a new graph is opened for every new measurement.

*\* This feature can be helpful for a certain type of experiment when the measured data are undesirably interpreted as negative values.*

---

## Icon Explanation:

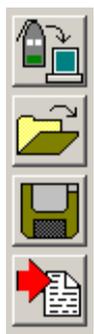

### Download

Downloads data from the AOM to PC.

### Load

Loads (opens) previously saved data files.

### Save

Saves data to hard disc.

### Export

Exports data in .txt format.

---

## 9.B. Example of Data Transfer and Visualization

**Step 1:** Perform a measurement with your AOM device (here, OJIP transient).

**Step 2:** Click the “Download” icon or select Device>Download.

**Step 3:** The Data table appears.

The screenshot shows the FluorPen software interface with a data table. The table has columns for Time, ID, and Measured parameter. The data is organized into three main sections, each with a callout:

- File name:** T:\Final\FluorPen\Test\_Data\data.dat
- Time of experiment:** 6:26:3 2.1 2007, 6:26:8 2.1 2007, 6:26:32 2.1 2007, 6:27:25 2.1 2007, 6:4:19 2.1 2007
- Measured parameter:** QY, OJIP, Ft, QY, OJIP

The table contains the following data:


Time	6:26:3 2.1 2007	6:26:8 2.1 2007	6:26:32 2.1 2007	6:27:25 2.1 2007	6:4:19 2.1 2007
ID	QY	OJIP	Ft	QY	OJIP
	<b>0.64</b>		<b>355</b>		<b>0.71</b>
Fo Backgr	1	Fo 618	Backgr 1	Fo Backgr 1	Fo 646
Fo Flash	351	Fj 1422	Flash 356	Fo Flash 275	Fj 1491
Fm Backgr	1	Fi 2149		Fm Backgr 1	Fi 2438
Fm Flash	964	Fm 2423		Fm Flash 942	Fm 2639
		Fv 1805			Fv 1993
		Vj 0.445			Vj 0.424
		Vi 0.848			Vi 0.899
		Fm/Fo 3.921			Fm/Fo 4.085
		Fv/Fo 2.921			Fv/Fo 3.085
		Fv/Fm 0.745			Fv/Fm 0.755
		Mo 0.911			Mo 0.831
		Area 4786890			Area 5229888
		Sm 2652.017			Sm 2624.129
		Ss 0.489			Ss 0.510
		N 5422.780			N 5142.671
		Phi_Po 0.745			Phi_Po 0.755
		Psi_o 0.555			Psi_o 0.576
		Phi_Eo 0.413			Phi_Eo 0.435
		Phi_Do 0.255			Phi_Do 0.245
		Phi_Pav 2857.311			Phi_Pav 2908.289
		ABS/RC 2.745			ABS/RC 2.595
		TRo/RC 2.045			TRo/RC 1.960
		ETo/RC 1.134			ETo/RC 1.129
		Dlo/RC 0.700			Dlo/RC 0.635
Description		Arabidopsis A			Arabidopsis B

Other callouts in the image include:

- Save experiment:** Points to the floppy disk icon in the left sidebar.
- Measured and calculated values:** Points to the 'Value' column in the table.
- Space for written notes:** Points to the 'Notes' tab at the bottom.
- Switch to graphic visualization of the experiment:** Points to the 'Graph:6:26:8 2.1 2007' tab at the bottom.
- Experiment description:** Points to the 'Description' column in the table.

**Step 4:** To visualize measurement in the graphic mode, click the “Graph” field in the bottom bar.

**Step 5:** The colored Graph of measured data appears.

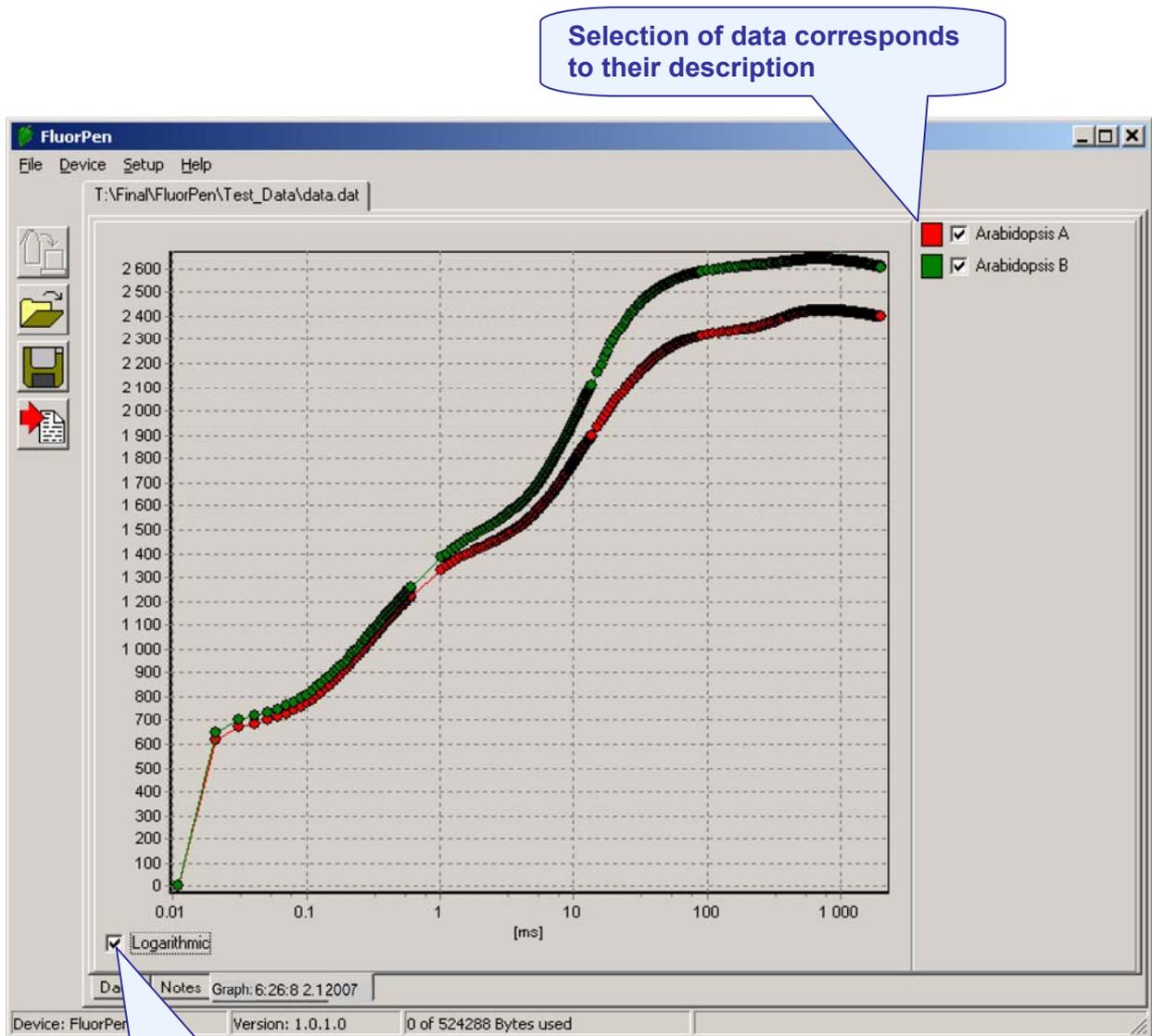

**Choose logarithmic or non-logarithmic graph form**

## 9.C. Explanation of OJIP Parameters

The AOM device offers the possibility to capture rapid fluorescence transient – OJIP, which occurs during exposure of photosynthesizing organisms to high irradiance. The FluorPen 2.0 software enables data downloading to a personal computer and subsequent OJIP analysis. The OJIP protocol includes the following extracted and technical parameters\*:

Formula Abbreviation	Formula Explanation
<b>Bckg</b>	<b>Background</b>
$F_0$	$F_0 = F_{50\mu s}$ , fluorescence intensity at 50 $\mu s$
$F_J$	$F_J$ = fluorescence intensity at J-step (at 2 ms)
$F_i$	$F_i$ = fluorescence intensity at i-step (at 60 ms)
$F_M$	$F_M$ = maximal fluorescence intensity
$F_V$	$F_V = F_M - F_0$ (maximal variable fluorescence)
$V_J$	$V_J = (F_J - F_0) / (F_M - F_0)$
$V_i$	$V_i = (F_i - F_0) / (F_M - F_0)$
$F_M / F_0$	
$F_V / F_0$	
$F_V / F_M$	
$M_0$ or $(dV/dt)_0$	$M_0 = TR_0 / RC - ET_0 / RC = 4 (F_{300} - F_0) / (F_M - F_0)$
<b>Area</b>	<b>Area between fluorescence curve and <math>F_M</math> (background subtracted)</b>
<b>Fix Area</b>	<b>Area below the fluorescence curve between <math>F_{40\mu s}</math> and <math>F_{1s}</math> (background subtracted)</b>
$S_M$	$S_M = \text{Area} / (F_M - F_0)$ (multiple turn-over)
$S_S$	$S_S$ = the smallest $S_M$ turn-over (single turn-over)
$N$	$N = S_M \cdot M_0 \cdot (1 / V_J)$ turn-over number $Q_A$
$\Phi_{P_0}$	$\Phi_{P_0} = 1 - (F_0 / F_M)$ (or $F_V / F_M$ )
$\Psi_{i_0}$	$\Psi_{i_0} = 1 - V_J$
$\Phi_{E_0}$	$\Phi_{E_0} = (1 - (F_0 / F_M)) \cdot \Psi_{i_0}$
$\Phi_{D_0}$	$\Phi_{D_0} = 1 - \Phi_{P_0} - (F_0 / F_M)$
$\Phi_{P_{av}}$	$\Phi_{P_{av}} = \Phi_{P_0} (S_M / t_{FM})$ $t_{FM}$ = time to reach $F_M$ (in ms)
$ABS / RC$	$ABS / RC = M_0 \cdot (1 / V_J) \cdot (1 / \Phi_{P_0})$
$TR_0 / RC$	$TR_0 / RC = M_0 \cdot (1 / V_J)$
$ET_0 / RC$	$ET_0 / RC = M_0 \cdot (1 / V_J) \cdot \Psi_{i_0}$
$DI_0 / RC$	$DI_0 / RC = (ABS / RC) - (TR_0 / RC)$

\* Formulas Derived From:

R.J. Strasser, A. Srivastava and M. Tsimilli-Michael (2000): The fluorescence transient as a tool to characterize and screen photosynthetic samples. In: *Probing Photosynthesis: Mechanism, Regulation and Adaptation* (M. Yunus, U. Pathre and P. Mohanty, eds.), Taylor and Francis, UK, Chapter 25, pp 445-483.

## 9.D. Software Update

---

---

### **Very important!**

**All data in the AOM memory are erased during the software update!**

**Before starting any software update, export all your data from the AOM memory into your computer!**

### Step 1: Starting Update


**Select:** Setup>Update Firmware



### Step 2: Selecting .bxn File

**Find:** Binary file (with the extension .bxn)

**Select:** Open

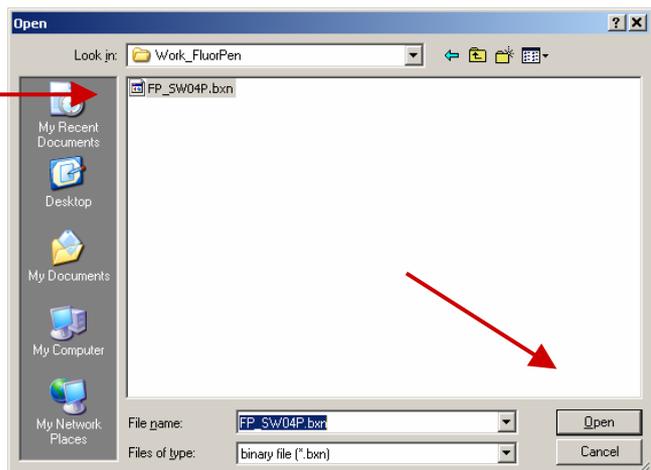

---

### Step 3: Finishing Upload

**Select:** “OK” to start uploading of the update.



The bottom bar indicates the upload progress.



**Press:** “OK” to finish upload.



## 10. Statement of Limited Warranty

---

- This Limited Warranty applies only to the AOM device and its accessories (excluding any batteries). It is valid one year from the date of shipment.
- If at any time within this warranty period the instrument does not function as warranted, return it and the manufacturer will repair or replace it at no charge. The customer is responsible for shipping and insurance charges (for the full product value) to PSI. The manufacturer is responsible for shipping and insurance on return of the instrument to the customer.
- No warranty will apply to any instrument that has been (i) modified, altered, or repaired by persons unauthorized by the manufacturer; (ii) subjected to misuse, negligence, or accident; (iii) connected, installed, adjusted, or used otherwise than in accordance with the instructions supplied by the manufacturer.
- The warranty is return-to-base only, and does not include on-site repair charges such as labor, travel, or other expenses associated with the repair or installation of replacement parts at the customer's site.
- The manufacturer repairs or replaces faulty instruments as quickly as possible; the maximum time is one month.
- The manufacturer will keep spare parts or their adequate substitutes for a period of at least five years.
- Returned instruments must be packaged sufficiently so as not to assume any transit damage. If damage is caused due to insufficient packaging, the instrument will be treated as an out-of-warranty repair and charged as such.
- PSI also offers out-of-warranty repairs. These are usually returned to the customer on a cash-on-delivery basis.
- *Wear & Tear Items* (such as sealing, tubing, padding, etc.) are excluded from this warranty. The term *Wear & Tear* denotes the damage that naturally and inevitably occurs as a result of normal use or aging even when an item is used competently and with care and proper maintenance.

*For customer support, please write to: [support@psi.cz](mailto:support@psi.cz)*

*Copyright © Photon Systems Instruments, 2009-10*